



FAQs for the stimulus preference chip

ScreenChip system compatibility

Do you need the ScreenChip system to operate the stimulus preference chip?

The Stimulus preference chip is a standalone chip. You not need a ScreenChip system in order to use it.

Set-up

What are the minimal requirements for set-up?

To operate the stimulus preference chip, you will need:

- (4) 10 mL fluid-reservoir syringes with one-way stopcock and 20G needle stubs
- (1) 5 mL syringe with 17G blunt needle stub
- (4) PE/6 polyethylene tubing
- (3) PE/9 polyethylene tubing
- (4) 1.0 mm OD SS tube
- (3) 1.5 mm OD SS tubes
- (1) stereomicroscope
- (2) vacuum sources
- Tape

See [Stimulus Preference Chip User Guide](#) for more information about the set-up.

How do I set-up the chip?

See [Stimulus Preference Chip User Guide](#)

Loading

How do you load the chip?

See [Stimulus Preference Chip User Guide](#)

How long does it take to load the chip?

It takes about 5 min to load the chip with testing fluids at the start of an experiment. It takes about the same amount of time to wash and load a worm.



Does it need to be taped to bench for loading?

We recommend viewing the chip on a stereomicroscope. Taping the chip to the microscope stage is necessary to keep it from moving as the stopcocks are adjusted, worms are inserted and removed

How many worms can you load at a time (is there a waiting area like for the SCs)?

For more consistent behavioral results from worm to worm, we recommend loading only one worm at a time into the chip. This is because there could be progressive changes in metabolic state as worms in the chip wait for their turn to perform.

How does preparation of the worm differ between the ScreenChip and the preference chip?

The main difference is that worms are washed in bulk for the ScreenChip, but individually for the Preference Chip.

How do you load the stimuli?

Stimuli are loaded into syringes the feed into the chip See Stimulus Preference Chip User Guide for further details.

Usage

What stage of worms does the stimulus preference chip accommodate?

The stimulus preference chip is designed to accommodate 1 day old adults.

What type of stimuli can the chip deliver?

The Stimulus Preference Chip can deliver any water soluble compound as a stimulus in any of the four inputs.

The chip can be operated in spatial or temporal modes:

- In the spatial mode, the solutions in the left and right channels are different. For example, the left and right solutions could be NaCl at different concentrations, yielding a step-wise concentration gradient. Or, the left and right solutions could contain the same concentration of NaCl and KCl, respectively, to measure the worm's preference for Na⁺ and K⁺ ions.
- In the temporal mode, the worm experiences a sudden shift between two stimuli that are each delivered symmetrically. For example, the left and right channels could first contain a solution low in NaCl that is then switched to a solution high in NaCl to measure the worm's response dynamics.

How many worms can I screen with one chip?



If solutions are well filtered to prevent the chip from clogging (e.g., 0.2 μm pore size), there is essentially no limit on the number of worms that can be tested in a single experimental session (e.g. one day in the lab). The chip can often be reused on subsequent days if it is cleaned and dried according standard procedures described in the [Stimulus Preference Chip User Guide](#). This procedure works best when inorganic compounds are used as stimuli. In the case of organic odorants, there is the concern that these are retained by the chip to an unknown degree after cleaning. In this case, it is safest to use a new chip between daily sessions. A similar concern exists when bacteria are used as stimulants.

Data

What type of data can be acquired with this chip?

The chip is designed primarily for the collection of behavioral data in several forms. These include the percentage of time the worm's head resides on the left or the right side of the chip, and mean head angle in the chip (measured with respect to the worm longitudinal axis of the body in the worm clamp).

How can I analyze the data obtained from the Stimulus Preference Chip?

Video recordings of worm behavior (30 frames/sec) can be analyzed in MATLAB using a custom routine to compute head angle θ in each image.

- 1- Frames mask and threshold frames to obtain an image of the worm.
- 2- Use a skeletonization procedure obtain the centerline of the worm.
- 3- Starting at the position of the restraint, traverse the centerline to find the tip of the head, defined as the point furthest from the restraint.
- 4- Initiation of reversals can be scored manually by an observer who was blind to experimental condition. Reversal behavior was defined as propagation of the undulatory wave in the posterior to anterior direction as previously described.

How long can I keep the worm "immobilized" in the chip?

You can keep the worm trapped in place for up to 20 min.

Can I recover my worms?

Yes, worms can be readily flushed out of the chip and recovered at the end of each experiment.